

Title:	Lifetime Turn Belt	ID:	
		0156	
Date in:	Response:	Model:	Author:
2005-12-07	2005-12-07	STX40 / STX200	CMa

Q:

I was asked by one of our Service Engineers to ask you about the shovel belts for the STX40 and STX200 incubators. He has seen where the current belt being used basically melts away in approximately 6 months in a system with Humidity and CO2. Is there another type of belt that we can use? He was under the impression that Cosmas stated that there was. Thanks in advance and have a great day.

A:

- 70°C shortens their lifetime.
- On the STX200 and on (very old) STX40s the motor torque could be set too high causing additional wear