

Title:	Handling does not react on Door Switch			ID:
				0372
Date in:	Response:	Model:	Author:	
2011-08-18	2011-08-18		CMA	

Q:

The current FW does not stop operation when glass door open we added special connector to these units so when the large enclosure door is open on the complete system the Freeslate switch would be connected to our units and trigger the same function as our glass door open , all should finish last operation and 1915 should be 0 ...

Customer called and told me that when glass door is open our units load unload plates anyway

Is there jumper or something that should have been removed / connected on this special connector in these units ? is it the FW ?

A:

Finishes operation - 1915 stays '0'. Door sw. 005 (1811) controls carrousel only. STR44 has no cars.

If you're suspicious about the hardware do

Check 1811 (005)

Check bypass toggle switch on module