

Title:	Difference Data to Display Value			ID:
				<i>0101</i>
Date in:	Response:	Model:	Author:	
<i>2010-09-21</i>	<i>2010-09-21</i>	-	<i>CMa</i>	

Q

The Display panel read 99.9% but the computer does not read above ~ 90.5% as seen below.

A

Could be broken display or broken N2 sensor. If sensor is broken display typically shows 99.9 but this value is also sent to controller.

Check display

1. Check if the display is still sending data. Open door and check if value on the PC changes.
2. Do RD DM995 (ADC N2). Overwrite this value WR DM995 9500 and read again RD DM995 to see if 9500 stays.