

Title:	Wiring Center Magnet of Tranferstation			ID:
				0257
Date in:	Response:	Model:	Author:	
2006-09-08	2006-09-08	-	CMa	

**Q:**

How is the Center Magnet on the Transfer Station wired?

**A:**

The Magnet is on Pin 3 on the 15-Pin xFer Strn. connector. On the 37-Pin connector is on Pin 32.

If you're sure they will never have a Swap Station you may, at the 37 Pin connector, move the wire on Pin 29 to Pin 32 and, at the 15 pin connector, move the wire from pin 9 to Pin 3.