

| | | | | |
|------------|--------------------------------|--------|---------|------|
| Title: | Sudden Temperature Drop | | | ID: |
| | | | | 0279 |
| Date in: | Response: | Model: | Author: | |
| 2006-10-16 | 2006-10-16 | - | CMa | |

Q:

If you look well at the file cooling 061006_2.bmp, you will see a drop of temperature at -23deg.C.

A:

By looking at the diagram it is hard to say what happened. I would first check the ventilation of the sensor. I added a pic. That shows the location of the sensor / vent. I also added a program STXMEX32 which allows you to check the performance of the chiller. Please reset the timers by pressing 'R' and confirm 'Y'. For the statistics the program does not need to run permanently. You may continue to use the system after the reset and then restart the program to read out the values 24h later.

