

Title:	Cassette Number of Positions			ID:
				0335
Date in:	Response:	Model:	Author:	
2012-12-13	2012-12-13	STX444	C <i>Ma</i>	

Q:

The STX44 should work with 2 cartages of 9 pos and two of 16 pos (the customer can change them), but the instrument can only work with the 9 pos , if we are using the 16 pos we get an error.

A:

The cassette pitch – which correlates with the plate height - is in DM23. The numbers of levels is in DM25 and must be updated as the number of levels change. The number of cassettes in DM9.