

<b>Title:</b>	<b>Defrost Temperature Peak</b>			<b>ID:</b>
				0364
<b>Date in:</b>	<b>Response:</b>	<b>Model:</b>	<b>Author:</b>	
2014-08-04	2014-08-04		CMa	

**Q:**

There is the peak between July 25 12:30 to July 26 14:00. We started to see such peak of the temperature in early-July > and it appears after defrost cycle. We didn't seen such peak last July. What do such peaks indicate? Is there anything we have to check?

**A:**

Yes the peak comes from a defrost cycle. Please log over longer period and also check how much water is coming on every defrost. If there is no or little water coming we may consider to increase cycle time.