

Title:	Lubrication Grease Z-Rail			ID:
				0378
Date in:	Response:	Model:	Author:	
2014-07-01	2004-04-27		C <i>Ma</i>	

Q:

Klübersynth LP 42-32 (for non-deep-freezing applications)

any other grease for rolling applications that is acid free works too. Pls. make sure to clean first with Washing Benzene (or petroleum ether). You may clean and grease the rack and the gears of the lift wagon too.

A:

Hi, my mate in Madrid is thinking to add some grease to Z guide in order to smooth the movement, do we can add grease? Let me know the right or recommended by Liconic, that can be also used in StoreX?